

Cellophane “A”, Cellulose, Plastics, Resins continued

Material	Moisture Range %	Autotransformer Setting	Infrared Lamp (w)	Drying Time (m)	Details
Resins, Paper, Fiber, Abrasives, Solid Glue	1.7-10.4	30-50 Depending on Material	250	10-15	
Resins, Plastic Powder	0.2-21.1	90-120	125	5-8	Composite report, samples finely ground
Resins, Vinyl	72	120	125	*	

Glauber Salt, Soap, Sulfanamides

Material	Moisture Range %	Autotransformer Setting	Infrared Lamp (w)	Drying Time (m)	Details
Glauber Salt – crude from filters	0-10	120	125	12-15	
Orthopedic Bandages	--	--	--	--	Not data provided, except use is satisfactory
Soap	10-14	120	125	10	
Soap Industrial Liquid	60-66	120	125	15	
Tablet Granulation type not specified	4-5	120	125	10	
Tablet Granulation of Triple Sulfanamides	0-30	120	125	7	
Terramycin – Culture in liquid media	95-99	120	125	15	